ENGLISH

NEED ASSISTANCE? Contact Celestron Technical Support
celestron.com/pages/technical-support

Product design and specifications are subject to change without prior notification. • This product is designed and intended for use by those 14 years of age and older.

CELESTRON ©2025 Celestron. Celestron and Symbol are trademarks of Celestron, LLC. • All rights reserved. • Celestron.com
US: Celestron, 2835 Columbia Street, Torrance, CA 90503 USA • UK: Celestron Global Ltd., Unit 2 Transigo, Gables Way, Thatcham RG19 4JZ, United Kingdom • Made in China | 05-25

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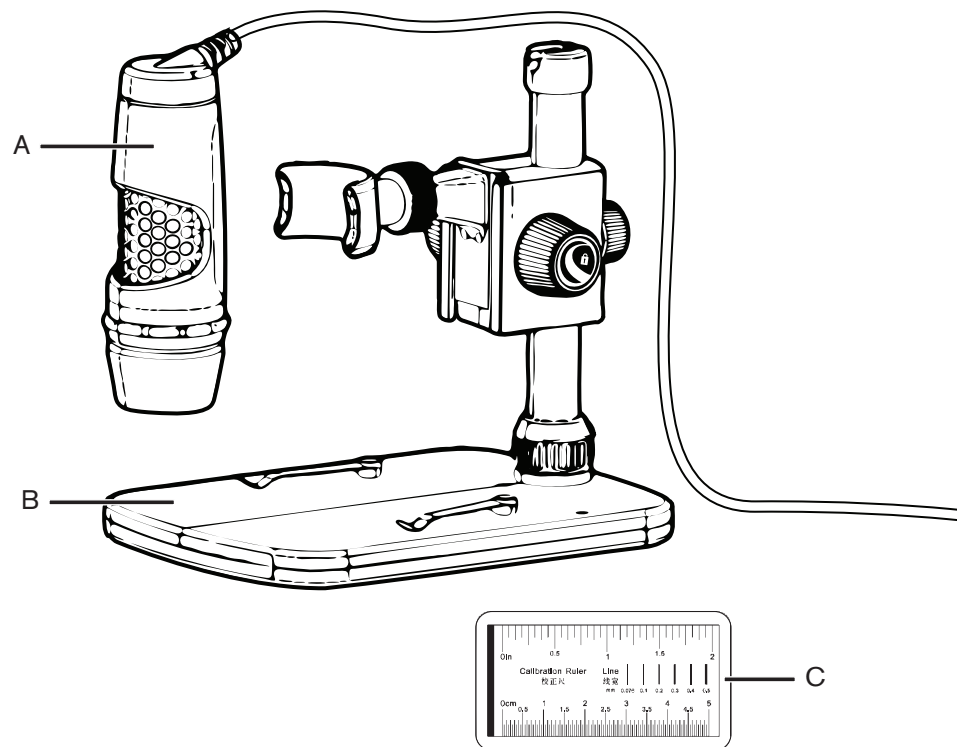
MODEL #44308

In the box

- [A] 1 x Celestron Labs 5MP USB-C digital microscope
- [B] 1 x Portable stand
- [C] 1 x Calibration ruler

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- [A] 1 x 5MP USB-C Digitalmikroskop
- [B] 1 x tragbarer Ständer
- [C] 1 x Kalibrierungslinéal

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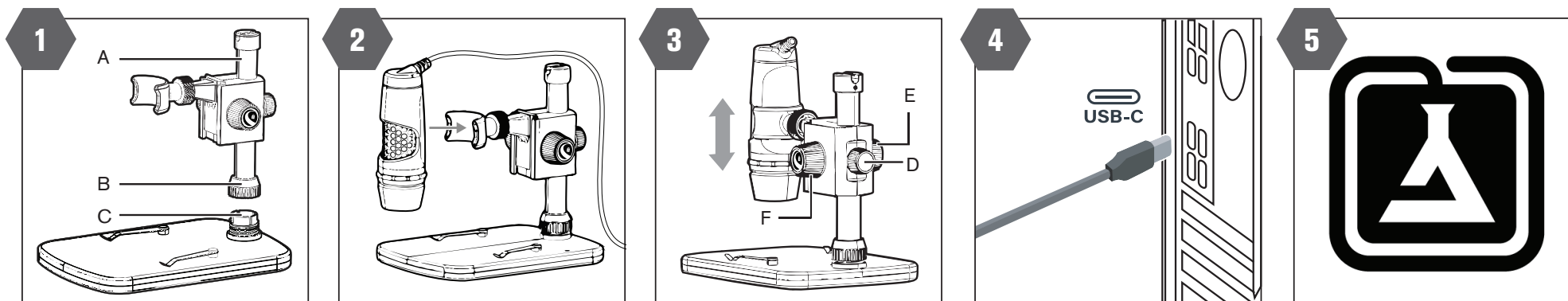
ENGLISH

1. Assemble the Stand

- Put the metal pole [A] through the coupler [B].
- Thread the coupler to fix the pole on the base [C].

2. Attach Microscope

- Snap the microscope into the U-shaped holder on stand.

3. Adjust Height

- Adjust the microscope's height by rotating main adjustment knob [D] and the side adjustment knob [E].
- Lock the microscope into place by rotating the lock knob [F].

4. Connect to Your PC

- Connect your handheld digital microscope via your computer's USB-C port.

5. Download Software & Manual

- On your PC, navigate to www.celestron.com/44308
- Download the Celestron Labs Digital Imaging Software.
- Download the full product manual.
- Launch the software to begin using your microscope.

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