For full manual visit [www.celestron.com](http://www.celestron.com)

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## ENGLISH

**NEED ASSISTANCE?** Contact Celestron Technical Support [celestron.com/pages/technical-support](http://celestron.com/pages/technical-support)

Product design and specifications are subject to change without prior notification. • This product is designed and intended for use by those 14 years of age and older.

**CELESTRON** ©2025 Celestron. Celestron and Symbol are trademarks of Celestron, LLC. • All rights reserved. • Celestron.com  
US: Celestron, 2835 Columbia Street, Torrance, CA 90503 USA • UK: Celestron Global Ltd., Unit 2 Transigo, Gables Way, Thatcham RG19 4JZ, United Kingdom • Made in China | 05-25

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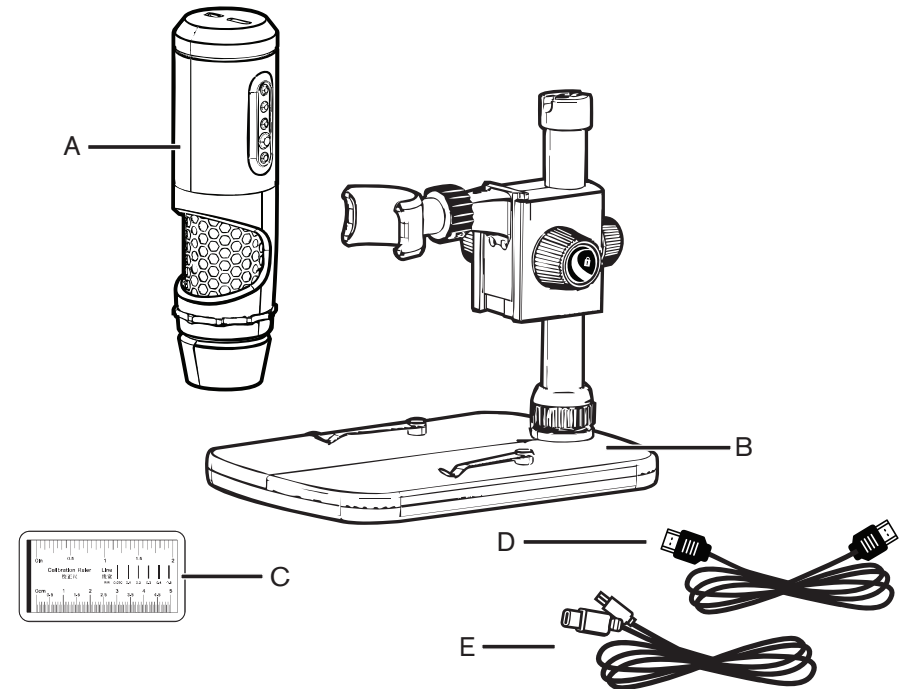
MODEL #44316

### In the box

- [A] 1 x MicroDirect® 1080p HD Microscope
- [B] 1 x Portable Stand
- [C] 1 x Calibration Ruler
- [D] 1 x HDMI Cable
- [E] 1 x USB-C Cable

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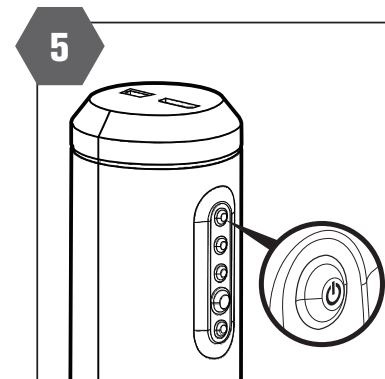
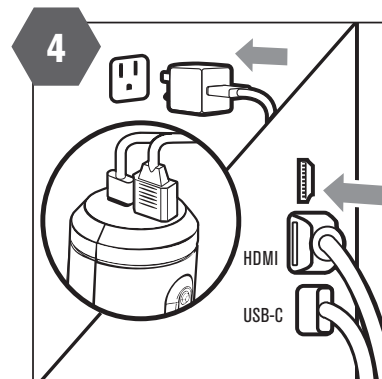
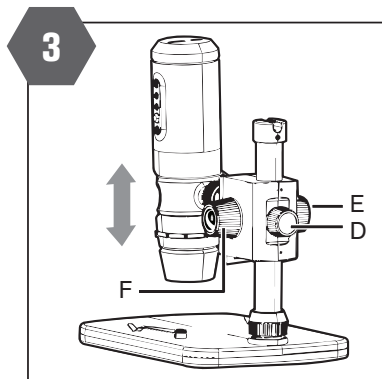
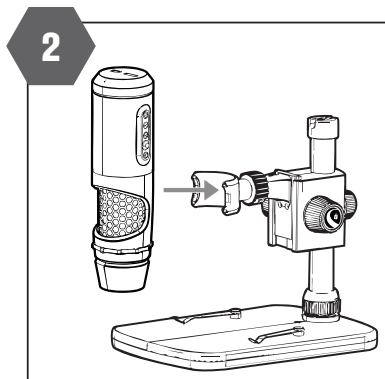
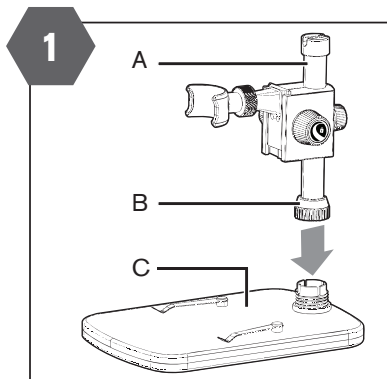
#### ENGLISH

### 1. Assemble the Stand

- Put the metal pole [A] through the coupler [B].
- Thread the coupler to fix the pole on the base [C].

### 2. Attach the Microscope

- Snap the microscope into the U-shaped holder on stand.

### 3. Adjust Height

- Adjust the microscope's height by rotating main adjustment knob [D] and the side adjustment knob [E].
- Lock down by rotating knob [F].

### 4. Connect Cables

- Connect one end of USB-C cable to microscope and the other end to a power 5V source.
- Connect one end of HDMI cable to microscope and the other end to a monitor, TV or projector.

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