**CLICK FOR FULL MANUAL**  
 CLIQUEZ POUR CONSULTER LE MODE D'EMPLOI COMPLET  
 FÜR DIE VOLLSTÄNDIGE ANLEITUNG KLICKEN  
 FARE CLIC PER IL MANUALE COMPLETO  
 PULSE PARA EL MANUAL COMPLETO

## ENGLISH

**NEED ASSISTANCE?** Contact Celestron Technical Support  
[celestron.com/pages/technical-support](http://celestron.com/pages/technical-support)

Product design and specifications are subject to change without prior notification. • This product is designed and intended for use by those 14 years of age and older.

Celestron hereby declares that the radio equipment type 44429 WiFi Micro Imager 2MP complies with Directive 2014/53/EU. The full text of the EU declaration of conformity is available at the following internet address [www.celestron.com/productcompliance](http://www.celestron.com/productcompliance).

**CELESTRON** ©2026 Celestron. Celestron and Symbol are trademarks of Celestron, LLC. • All rights reserved. • Celestron.com  
 US: Celestron, 2835 Columbia Street, Torrance, CA 90503 USA • UK: Celestron Global Ltd., Unit 2 Transigo, Gables Way, Thatcham RG19 4JZ, United Kingdom • EU Authorized Representative: AR Experts B.V., Boeingavenue 209, 1119 PD Schiphol-Rijk, The Netherlands, info@ar-experts.eu • Made in China • 03-26

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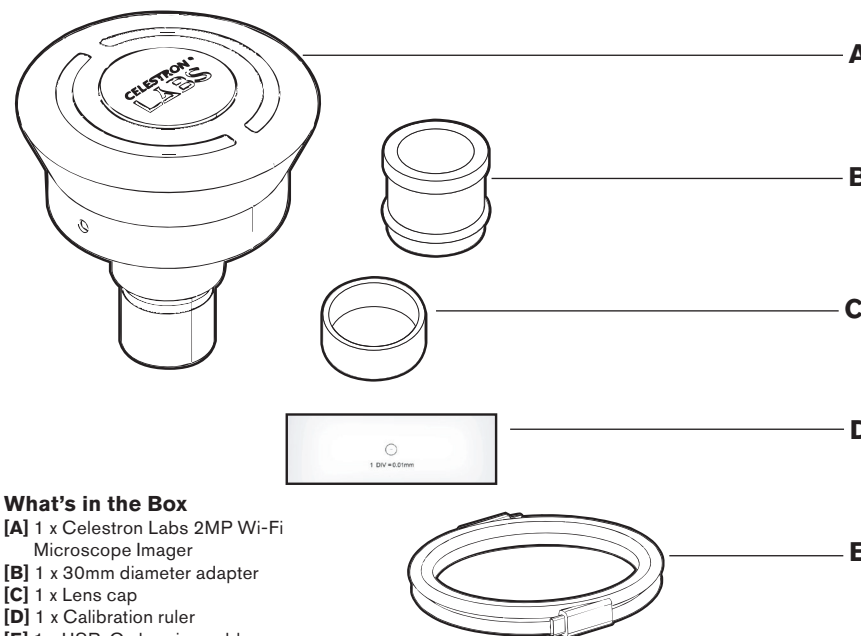
## QUICK START GUIDE

### What's in the Box

- [A] 1 x Celestron Labs 2MP Wi-Fi Microscope Imager
- [B] 1 x 30mm diameter adapter
- [C] 1 x Lens cap
- [D] 1 x Calibration ruler
- [E] 1 x USB-C charging cable

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
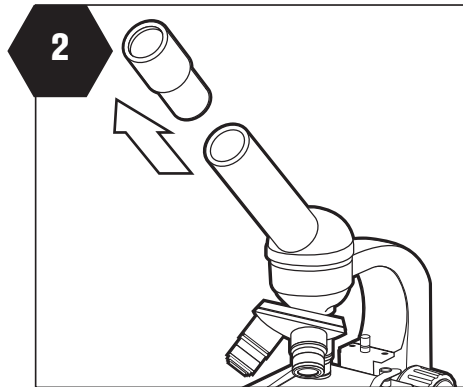
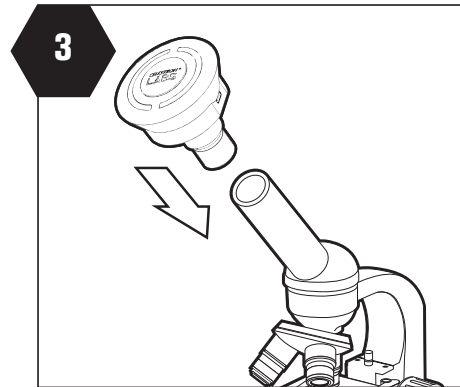
**3. Place the imager into the microscope barrel.**

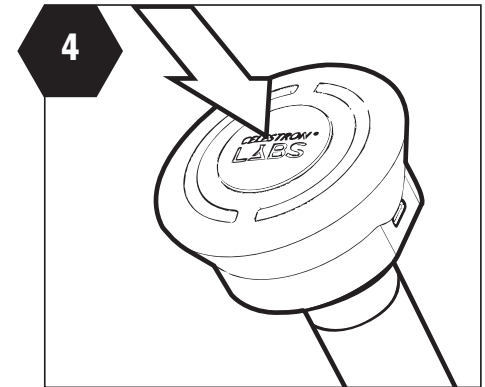
- 3. Placez l'imageur dans le corps du microscope.
- 3. Setzen Sie den Imager in das Mikroskopgehäuse ein.
- 3. Inserire l'imaginer nel barilotto del microscopio.
- 3. Coloque el generador de imágenes en el cuerpo del microscopio.

**4. Press the power button to turn on the imager.**

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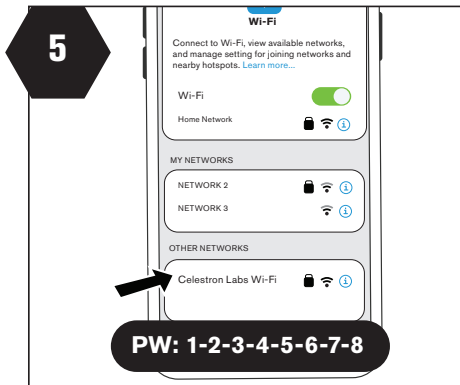
**5. Connect with Celestron Wi-Fi network through your smart device.**

- 5. Connectez-vous au réseau Wi-Fi Celestron via votre appareil intelligent.
- 5. Stellen Sie über Ihr Smart-Gerät eine Verbindung mit dem Celestron-WLAN-Netzwer
- 5. Connettiti alla rete Wi-Fi Celestron tramite il tuo dispositivo smart.
- 5. Conéctese a la red Wi-Fi de Celestron a través de su dispositivo inteligente.


**6. Open the Celestron Labs Wi-Fi app.**

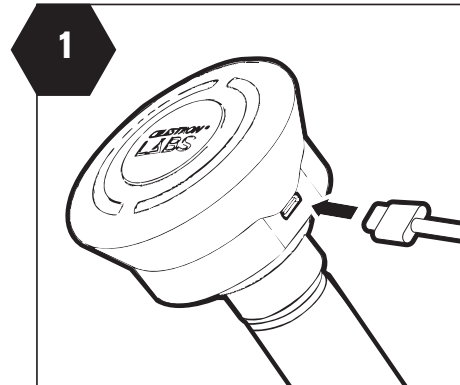- 6. Ouvrez l'application Wi-Fi Celestron Labs.
- 6. Öffnen Sie die Celestron Labs Wi-Fi-App.
- 6. Aprire l'app Wi-Fi di Celestron Labs.
- 6. Abra la aplicación Wi-Fi de Celestron Labs.

**1. Connect one end of USB-C cable (E) to the imager.**

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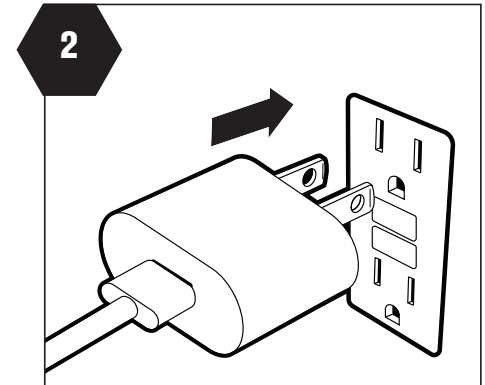**2. Connect one end of USB-C cable (E) to your 5V-1.0 A to 1.8A AC adapter and plug in.**

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